



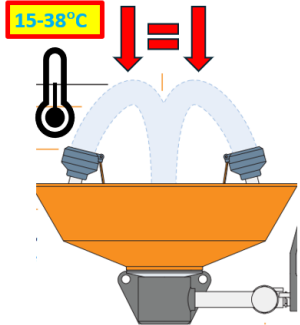




# Emergency Eyewash Testing

1		<ul style="list-style-type: none"> <li>➤ Test eyewash <b>weekly</b>.</li> </ul>
2		<ul style="list-style-type: none"> <li>➤ Check that area around eyewash is <b>unobstructed and uncluttered</b>.</li> </ul>
3		<ul style="list-style-type: none"> <li>➤ <b>Activate</b> the unit.             <ul style="list-style-type: none"> <li>○ Place a shield or a 5-gallon bucket under it, if needed.</li> </ul> </li> </ul>
4		<ul style="list-style-type: none"> <li>➤ Check that protective eyewash <b>covers come off</b> readily.</li> </ul>
5		<ul style="list-style-type: none"> <li>➤ Check for <b>constant, even flow</b> from both eyewash pieces, <b>appropriate pressure</b> (to reach eyes), and <b>tepid temperature</b> (15-38°C).</li> </ul>
6		<ul style="list-style-type: none"> <li>➤ <b>Run water for a few minutes</b> and check that the drain works.</li> </ul>
7		<ul style="list-style-type: none"> <li>➤ Mark the <b>inspection tag</b>.</li> <li>➤ <b>Report any problem</b> to Facilities immediately to have it corrected.</li> </ul>